

Version No.: FAL1715

GenFQ Hot Start Taq Master Mix

GenFine Code: A109-01, A109-02

Storage condition: Transportation at $\leq 0^{\circ}\text{C}$, Store at $-30 \sim -15^{\circ}\text{C}$, Valid period 2 years.

Description:

GenFQ Hot Start Taq Master Mix is a $2 \times$ concentrated solution, including GenFQ Hot Start Taq DNA polymerase, dNTPs and all other components required for PCR, except DNA template and primers. This pre-mixed formulation saves time and reduces contamination due to a reduced number of pipetting steps required for PCR set up. The mix is optimized for efficient and reproducible PCR. The protective agent added in the amplification system made $2 \times$ Taq Master Mix remained stable activity after repeated freezing and thawing. It has the advantages of high sensitivity, specificity and stability.

Components:

Components	A109-01	A109-02
$2 \times$ Hot Start Taq Master Mix	1 ml	1 ml $\times 5$
ddH ₂ O	1 ml	1 ml $\times 5$

Applications:

Gene detection: especially suitable for large-scale gene detection, semi quantitative PCR experiment and micro DNA detection, High specificity, sensitivity PCR reaction, Genomic amplification with strong background (such as the detection of a specific gene site or foreign pathogen in the genome), etc.

Protocol:

1. Prepare the reaction solution as follows:

Components	Volume (50 μl)
$2 \times$ Hot start Taq Master Mix	25 μl
Forward Primer (10 μM)	1 μl
Reverse Primer (10 μM)	1 μl
Template DNA*	X μl
ddH ₂ O	Up to 50 μl

* Plants and animals Genomic DNA 1~500 ng, E.coli Genomic DNA 1~100 ng, λ DNA 0.1~10 ng, Plasmid 0.1~10 ng.

2. Thermal cycling conditions:


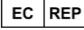












Perform PCR using recommended thermal cycling conditions:

Step	Temperature	Time	Number of cycles
1	95 $^{\circ}\text{C}$	2 min	1 cycle

2	95°C	15~30 sec	25~35 cycles
	55°C*	15~30 sec	
	72°C	1 kb/min	
3	72°C	5~10 min	1 cycle

*The annealing temperature should be 3~5°C lower than the melting temperature (T_m) of the primers. For complex template, the annealing temperature must also be adjusted.

[Symbols]

Symbols	Meanings
	Manufacturer
	Authorized representative in the European Community
	<i>In vitro</i> diagnostic medical device
	This product fulfills the requirements of the European Directive 98/79 EC for <i>in vitro</i> diagnostic medical devices.
	Catalogue number
	Batch code
	Date of manufacture
	Use-by date
	Temperature limit
	Consult instructions for use
	Keep dry
	Keep away from sunlight
	Do not re-use
	Do not use if package is damaged



GENFINE BIOTECH (CHANGZHOU) CO., LTD.

4th Floor, Building E4, No.9, Changyang Road, West Taihu Technology Industrial Park, Changzhou City, Jiangsu Province

PEOPLE'S REPUBLIC OF CHINA

Tel: +86 051983761557

E-mail: marketing@genfine.com

Web: en.genfine.com



Lotus NL B.V.

Koningin Julianaplein 10, 1e Verd, 2595AA, The Hague, Netherlands.

E-mail: peter@lotusnl.com

Tel: +31644168999