

Version No.: FDL1715

GenFQ Taq DNA Polymerase

GenFine Code: A111-01, A111-02

Storage condition: Transportation at $\leq 0^{\circ}\text{C}$, Store at $-30 \sim -15^{\circ}\text{C}$, Valid period 2 years.**Description:**

GenFQ Taq DNA Polymerase is a highly thermostable recombination DNA polymerase with a $5' \rightarrow 3'$ DNA polymerase activity and a $5' \rightarrow 3'$ exonuclease activity, and lacks a $3' \rightarrow 5'$ proofreading function. This product can be used in a variety of templates and has good tolerance to fluorescent labeled dUTP and dITP nucleotides. The 3'-end of the PCR products contain A, which can be directly cloned to a T-vector.

Components:

Components	A111-01 (500 U)	A111-02 (1000 U)
GenFQ Taq DNA Polymerase	100 μl (5 U/ μl)	200 μl (5 U/ μl)
5 \times Taq Buffer	1.1 ml	1.1 ml*2

Applications:

Widely used for DNA amplification of animals, plants and microorganisms.

Definition of Activity Unit:

One unit of the enzyme catalyzes the incorporation of 10 nmol of deoxyribonucleotides into a polynucleotide fraction in 30 min at 74°C .

Certificate of Analysis:

- 1) Endodeoxyribonuclease Assay: No conversion of covalently closed circular DNA to nicked DNA was detected after incubation of 10 U of Taq DNA Polymerase with 1 μg of pUC19 DNA for 4 hours at 37°C .
- 2) Exodeoxyribonuclease Assay: No degradation of DNA was observed after incubation of 1 μg of lambda DNA/Hind III fragments with 10 U Taq DNA Polymerase for 4 hours at 37°C .
- 3) Detection of E. coli DNA residue: 20 μl the residual nucleic acid in the product was detected by TaqMan qPCR method, and the residue of E. coli genome was less than 10 copies.

Protocol:**1. Prepare the reaction solution as follows:**

Components	Volume (50 μl)
Template DNA*	X μl
GenFQ Taq DNA Polymerase	1 μl
Forward Primer (10 μM)	1 μl
Reverse Primer (10 μM)	1 μl
5 \times Taq Buffer	10 μl
dNTPs (10 mM)	1 μl
ddH ₂ O	Up to 50 μL

* Plants and animals Genomic DNA 1~500 ng, E.coli Genomic DNA 1~100 ng, λ DNA 0.1~10 ng, Plasmid 0.1~10 ng.

2. Thermal cycling conditions:

For research use only, not for use in diagnostic procedures.

Perform PCR using recommended thermal cycling conditions:


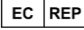












Step	Temperature	Time	Number of cycles
1	95°C	2 min	1 cycle
2	95°C	15~30 sec	25~35 cycles
	55°C*	15~30 sec	
	72°C	1 kb/min	
3	72°C	5~10 min	1 cycle

*The annealing temperature should be 3~5°C lower than the melting temperature (T_m) of the primers. For complex template, the annealing temperature must also be adjusted.

Notes:

Taq DNA Polymerase has a certain reaction activity at room temperature, the PCR reaction system should be prepared on ice, and then placed on the PCR instrument for reaction. This can reduce the non-specific amplification in the reaction preparation stage and help to obtain high specific amplification results.

[Symbols]

Symbols	Meanings
	Manufacturer
	Authorized representative in the European Community
	<i>In vitro</i> diagnostic medical device
	This product fulfills the requirements of the European Directive 98/79 EC for <i>in vitro</i> diagnostic medical devices.
	Catalogue number
	Batch code
	Date of manufacture
	Use-by date
	Temperature limit
	Consult instructions for use
	Keep dry
	Keep away from sunlight
	Do not re-use
	Do not use if package is damaged



GENFINE BIOTECH (CHANGZHOU) CO., LTD.
 4th Floor, Building E4, No.9, Changyang Road, West Taihu Technology Industrial
 Park, Changzhou City, Jiangsu Province
 PEOPLE'S REPUBLIC OF CHINA
 Tel: +86 051983761557
 E-mail: marketing@genfine.com
 Web: en.genfine.com



Lotus NL B.V.
 Koningin Julianaplein 10, 1e Verd, 2595AA, The Hague, Netherlands.
 E-mail: peter@lotusnl.com
 Tel: +31644168999